




14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ [www.coylab.com](http://www.coylab.com)

# Compact Dehumidifier Instructions

## Dehumidifying module:

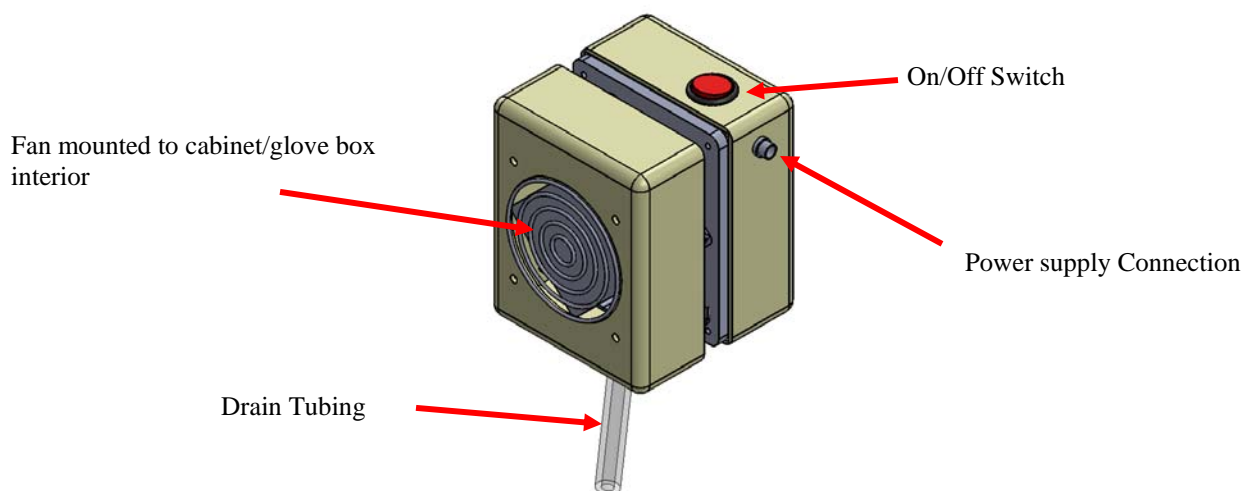

## How to use:

1. Turn on the dehumidifier (red led switch on outside top of the unit).
2. The unit is always on . A continuous forced circulation design ensures accurate and uniform humidity level inside the cabinet or glove box.
3. You will need to empty the water reservoir from time to time depending on how much water is collected.
4. To ensure efficient dehumidification, do not block the air circulation fans.

### Specifications:

- **Input Power:** DC 12V  $\pm$  10% / 6.67dc A / 60W
- **Humidity Range:** +25% Rh
- **Power supply:** AC 120-220 V 60Hz
- **1.224 pounds/0.551kg of water per 24 hours**