



14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ www.coylab.com

Calibrator retrofit sensor port for S15 02 sensor

Old Fitting:



1. using Channel locks you can remove the old sensor port fitting. (Counter Clockwise)





14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ www.coylab.com

2. Old fitting removed.



3. Shown with new larger fitting for the S15 O2 sensor.





14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ www.coylab.com

4. Screw larger fitting on and finger tight is good. Do not over tighten. An air tight seal is not required.



5. Below is how the S15 O2 sensor should be position when putting into the calibrator.

