



14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ www.coylab.com

Sensor Port Retrofit for Cabinet or Glove box

If you purchased a new O₂ sensor for hypoxic chamber/cabinets and the glove box/cabinet is 2 years or older you may require a new sensor adaptor in order to fit the new larger O₂ sensor. (pictured page 2) Please contact Coy at techservice@coylab.com or call 734-457-2200 and we can send you one.

1. Remove old fitting with crescent wrench or channel locks. (Counter clockwise)



2. Place new sensor port into hole with the O-ring on the bottom side of the sensor port.





14500 Coy Drive ▪ Grass Lake, MI 49240 ▪ Ph: (734) 475-2200 ▪ Fx: (734) 475-1846 ▪ www.coylab.com

3. using channel locks push down on the sensor port as you tighten with the locks. (clockwise)



4. Silicone stopper with predrilled hole and sensor in final location. Now sensor is inside the chamber at uniform temperature.

