

Meet the Coy Quick Change Cuff

Faster swaps. Fewer leaks. Less gas wasted.

A few years ago, we noticed a consistent user problem with our Coy chambers. Despite the importance of changing gloves once they began to show signs of rare, users would consistently avoid this task:

- The process could take up to half an hour.
- Replacement required attention and training.
- Too many users put off the task or left it to someone else.

However, delaying glove replacement can lead to increased gas consumption, higher operational costs, and a compromise of the anaerobic state within the chamber.



Coy's Quick Change Cuff System Reduces Swap Time to Under a Minute

The Coy Quick Change Cuff takes less than a minute to replace a glove with a simple twist and catch motion. Users seat the glove on the cuff, close off the cuff attachment from outside the chamber, twist off the existing cuff and twist on the new unit.

Since 2019 the Quick Change Cuff is standard on all new Coy Vinyl Anaerobic Chambers.

Older chambers can be retrofitted, however doing so may not be cost effective if the current chamber is nearing the end of its lifecycle.